


## Bacterial analysis in flowcytometry

Written by: Anja Bohn

Date: May 2020, modified Nov 2023

### Sample preparation

- To be able to differentiate between bacteria and background we highly recommend using a DNA stain
- It is very important to standardize the staining of your samples. Both incubation temperature and incubation time will affect the outcome. See figure 1.

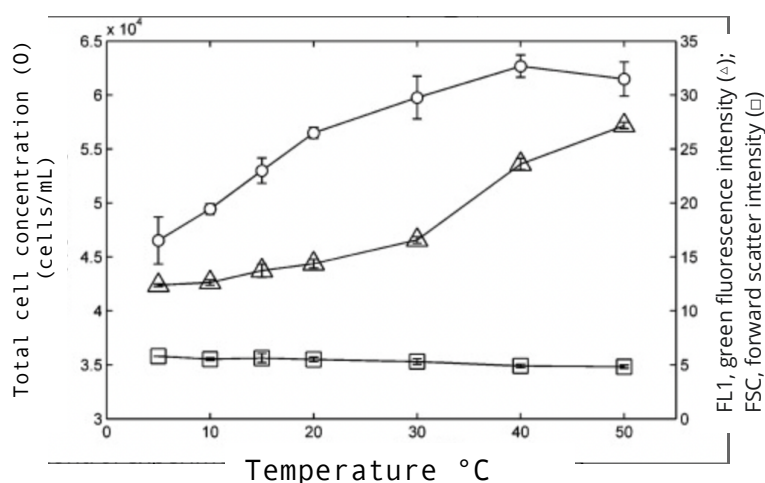


**Figure 1:** Effect of different incubation times and incubation temperature on bacteria count  
*Frederik Hammes, <https://onlinelibrary.wiley.com/doi/10.1002/cyto.a.22048>*

### Analysis

- The Quanteon and the 3-laser FACS Aria III are the best instruments for bacteria analysis and sorting in the FACS Core Facility.
- Use height (not area) for fluorescence as well as scatter signals. Because of the small size of bacteria, there is less variation on the height parameter compared to area.
- Use log scales on both forward and side scatter.
- If sorting on the FACS Aria III remember the possibility to remove the ND filter.
- Set your threshold on a fluorescence parameter (DNA stain). If this is not possible, a side scatter threshold is often better than a forward scatter threshold.

- Make sure you look at single events
  - Start out using the slowest speed possible (not just “slow”, but move the slider to the far left)
  - Determine the median fluorescence intensity (MFI) for the bacteria population.
  - Repeat this step but on a diluted version of your first sample.
  - The count-rate should decrease, and the MFI should stay the same as in the undiluted sample. If this is not the case, you are exposed to swarming (many events are being detected as one event). In this case you should make a dilution series to determine the dilution where you are detecting single events. See figure 2.

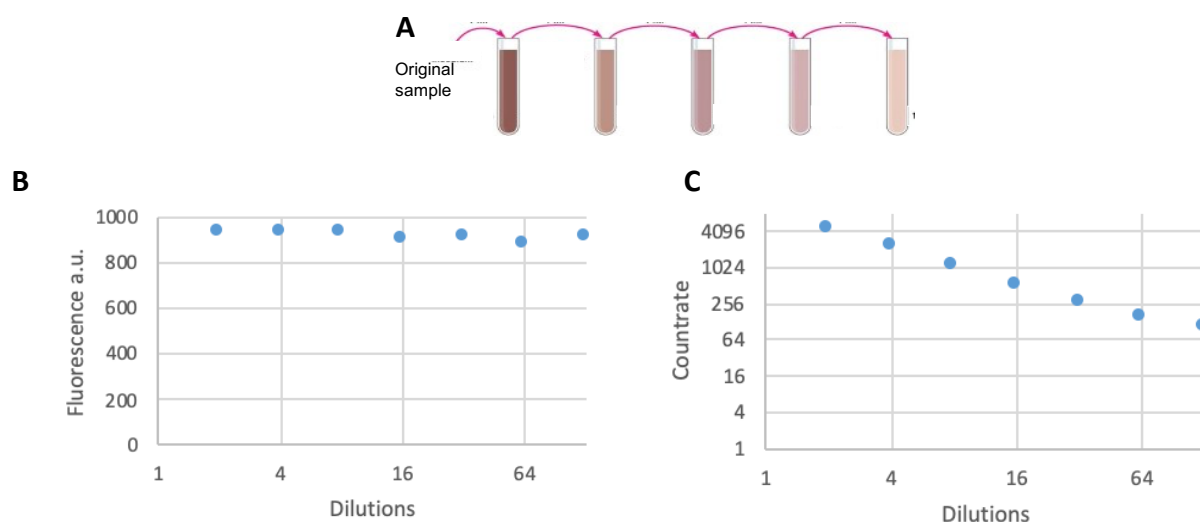

**Figure 2:** Dilution series of bacterial sample (A) does not affect fluorescence intensity (B), but the event-rate/count-rate decreases (C).

*Ref: Denis Selnihhin, Department of Molecular Biology and Genetics - Geneexpression and Genomedicine, Aarhus University and Anja Bille Bohn, FACS Core Facility*